AD-8129TH

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
Simplified Instruction Manual

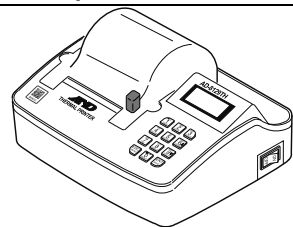View the product page

Refer to the detailed instruction manual on the product page: <https://link.aandd.jp/ad8129>

About this manual

- (1) No parts of this manual may be reproduced without permission. This manual may not be copied, modified, or translated without the written permission of A&D Company, Limited.
- (2) The contents of this manual are subject to change without notice.
- (3) Please contact A&D if you notice any uncertainty, errors, omissions, etc. in this manual.
- (4) A&D Company, Limited bears no liability for direct, indirect, special, or consequential damages due to the operation of the product, even if advised of the possibility of such damage. Furthermore, A&D assumes no liability for claims of rights from third parties. Concurrently, A&D assumes no liability whatsoever for data losses regardless of (3) above.

© 2024 A&D Company, Limited

A&D Company, Ltd.

1WMPD4005386


1. Introduction

This manual is a product overview of the AD-8129TH and instructions for assembly and printing. For details on operation and communication methods, please check the information on the A&D website.

2. Features

- This device is a printer used to connect to a weighing device such as A&D's electronic balance, scale, counting scale, moisture analyzer, vibro viscometer or weighing indicator and print out data transmitted from the weighing device.
- With thermal printing, the printed contents can be clearly and easily viewed, and the print density can also be changed.
- Light weight and compact size of the printer enables it to be set up in various locations and be easily transported.
- The printer is equipped with the RS-232C or current loop interface for data input, so it allows easy connections between the printer and a weighing device such as an electronic balance.
- Built-in clock function of the printer enables printing out of the date and time.
- Printing at a regular interval is possible. (Interval printing)
- Statistical calculation results of the weighing data, including standard deviation, can be printed out.
- Time series variation of the weighing data in chart form can be printed out.
- The liquid crystal display provides easy conditions confirmation and settings of the printer.
- By pressing the key switch of the printer, a re-zero command to zero the display and request commands for the weighing data can be transmitted to a weighing device, enabling the weighing data to be printed out without touching the highly sensitive weighing device.

3. Package contents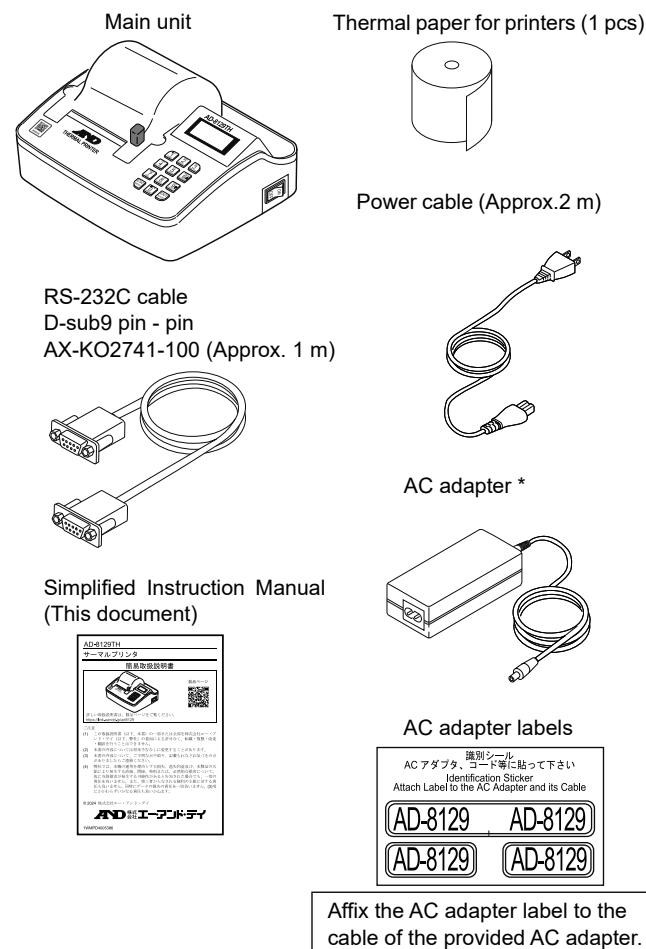


* Confirm that the AC adapter and power cable are correct for the local voltage and power receptacle type.

4. Assembly


□ Connecting the AC adapter

1. Confirm that the power switch is turned off.
2. Connect the power cable and AC adapter.
3. Connect the DC plug of the AC adapter to the power connector of this unit.
4. Connect the AC plug to the outlet.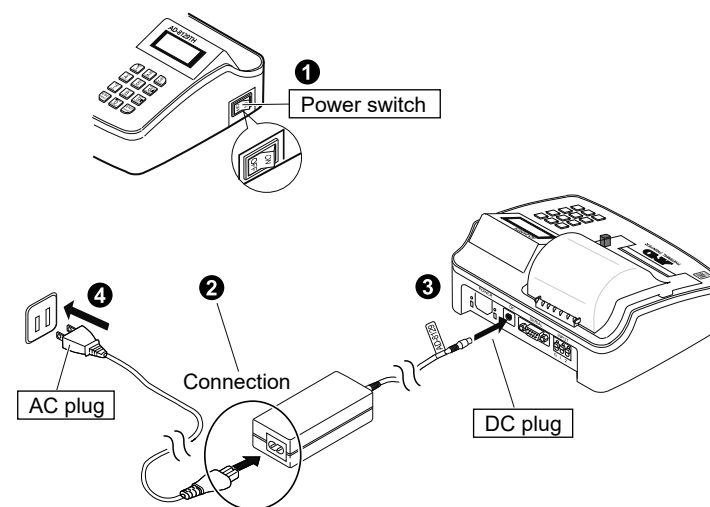


□ Opening and closing the printer cover

1. Press the button on the right side of the printer cover.
2. Lift the printer cover upwards.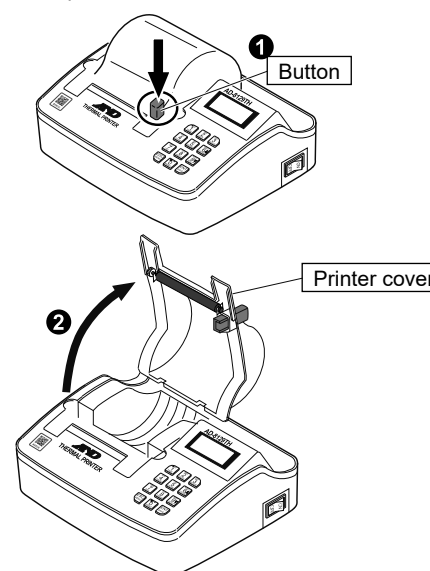

□ Installing the printer paper

1. Place the printer paper so that its leading edge protrudes from the printing unit.
2. Close the printer cover.
3. Turn the power switch on.
4. Press the key to feed the printer paper 5 or 6 cm out of the printing unit.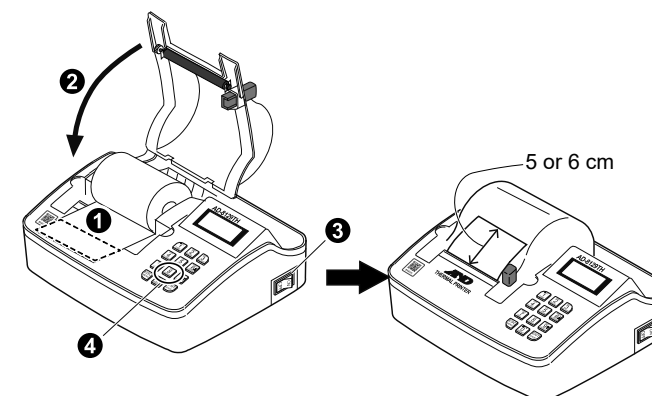


Caution

The AD-8129TH is a thermal printer designed exclusively for use with thermal paper. The plain paper for AD-8127 cannot be used.

□ Connecting to the weighing device

Use the RS-232C cable or current loop cable to connect the printer to the weighing device. The factory default communication settings are baud rate: 2400 bps, data length: 7 bits, parity: EVEN, stop bit: 1 bit. This is in accordance with the factory default settings for most of A&D's weighing devices.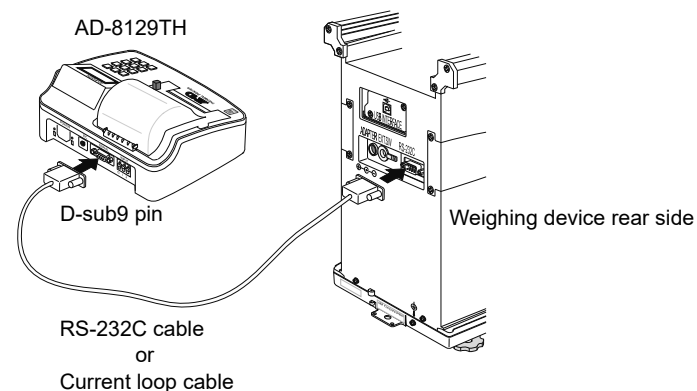

5. Printing

□ Printing out data from the weighing device

1. Turn the printer power on. The factory default printing mode is "External key printing mode (EXT. KEY)".
2. Connect the printer to the weighing device using a RS-232C or current loop cable.
3. Transmit data to the printer from the weighing device. (e.g.) Press the key on the weighing device.
4. Press and hold the key to feed the printer paper until the printed results can be viewed, and then cut off the printer paper.

□ Paper feed

Press the key to allow paper feed while the button is pressed.

For detailed instructions on how to use, please refer to the instruction manual on the product page.

6. Storage of thermal paper

- Store the thermal paper in a cool, dry, and dark place.
- Do not rub the thermal paper with hard objects.
- Do not leave the thermal paper near organic solvents.
- Do not allow the product to come into contact with PVC film, erasers, or adhesive tape to touch the thermal paper for extended periods of time.
- Do not stack the thermal paper with whiteprints or wet copies immediately after copying.

7. Specifications

Printing method	Thermal type
Character specifications	13 dots × 28 dots 1.5(W) × 2.6(H) mm
Number of characters printed	24 characters per line
Operating temperature range	0 °C to 40 °C
Operating humidity range	20 % to 85 %RH (No condensation)
Storage temperature range	-10 °C to 50 °C
Storage humidity range	10 % to 90 %RH (No condensation)
Interface	RS-232C, current loop
Power supply & AC adapter type	Confirm that the AC adapter and power cable are correct for the local voltage and power receptacle type.
External dimensions	174(W)×141(D)×87(H) mm
Weight	Approx. 500 g
Accessories	Thermal paper for printers (1 roll), Power cable, AC adapter, AC adapter labels, RS-232C cable (D-sub9 pin - 9 pin, 1 m), Simplified Instruction Manual

8. Sold separately

Thermal paper for printers (5 pcs)	AX-PP147-S
Dust-free thermal paper for printers (10 pcs)	AX-PP183-S
RS-232C cable (D-sub9 pin - 9 pin, 1 m)	AX-KO2741-100
RS-232C cable (D-sub9 pin - 9 pin, 2 m)	AX-KO2466-200
RS-232C cable (D-sub25 pin - 9 pin, 0.9 m)	AX-KO2116-090
RS-232C cable (D-sub25 pin - 9 pin, 2 m)	AX-KO1710-200
RS-232C cable (DIN - 9 pin, 2 m)	AX-KO1786-200
Current loop cable (DIN - 9 pin, 1 m)	AX-KO5291-100
Foot switch	AX-SW138