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
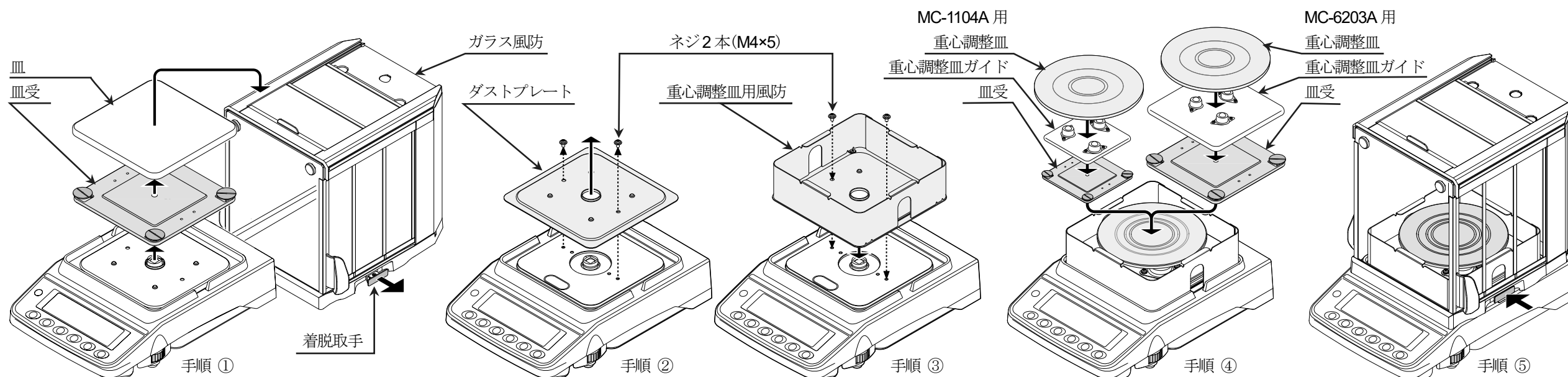
Auto-Centering Pan (AX-MC1104A-PAN / AX-MC6203A-PAN) Instruction Manual

Read this manual thoroughly before using the auto-centering pan (AX-MC1104A-PAN/AX-MC6203A-PAN) and keep it at hand for future reference.

1. Introduction

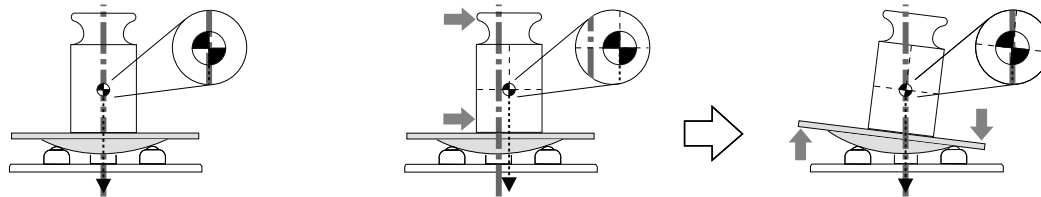The auto-centering pan can align the center of gravity of the material being weighed (a weight) with the pan center automatically and reduce eccentric loading errors when this pan is installed onto a MC-1104A or MC-6203A balance. The influence of drafts can be reduced by attaching the breeze break provided with this option.

Center of gravity of the weight and the pan are aligned.

When the center of gravity of the weight and the pan are not aligned, the center of gravity of the weight shifts to the center of auto-centering pan the automatically.

3. Installation

Step 1 Pull both **breeze break locking handles** to remove the **glass breeze break unit**.
Remove the **pan support** and **weighing pan**.

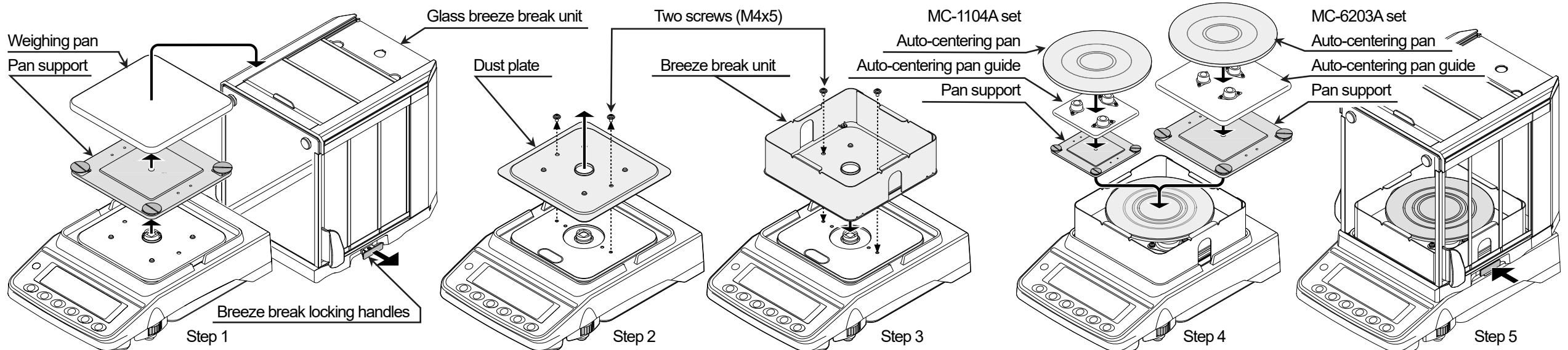
Step 2 Remove the **two screws** (M4x5) that secure the **dust plate** and remove the **dust plate** from the balance.

Step 3 Install the **breeze break unit** using the **two screws** (M4x5) removed at step 2.


Step 4 Place the **pan support** on the balance, place the **auto-centering pan guide** on the pan support, and then place the **auto-centering pan** on that.
While tilting (forward, backward, left and right) the **auto-centering pan guide** of the MC-6203A, put it on the **pan support**.

Step 5 To install the **glass breeze break unit**, place it on the balance and push in both **breeze break locking handles**.


CAUTION When replacing the weighing pan with the auto-centering pan, be sure to adjust sensitivity of the balance before weighing because the weights of the two pans are different.

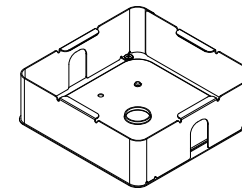2. Unpacking


Breeze break unit

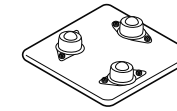
Auto-centering pan guide for MC-1104

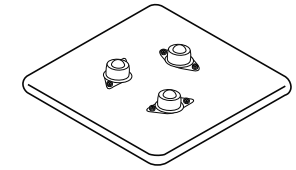Either pan guide is included in the packing.

Auto-centering pan guide for MC-6203

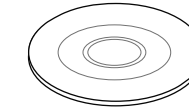Auto-centering pan

4. Precautions to use

- Do not place the material being weighed on the edge of the auto-centering pan. We recommend placing the material being weighed on the auto-centering pan in a manner where the pan is as horizontal as possible.
The auto-centering pan may touch the breeze break and cause the material being weighed to fall, or the center of gravity of the material being weighed may not shift to the pan center, causing a weighing error.
- Place a weight on the auto-centering pan gently.
Dropping a weight onto the auto-centering pan may damage to the surface of the pan where it contacts the auto-centering pan guide bearings, affecting smooth centering of the pan.

A&D
A&D Company, Limited