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
1. Features

The FX-11 protects the balance from the effects of a breeze and will provide more accurate weighing results. All components are antistatic, making it less likely that weighing errors will occur due to static electricity. When used with the FZ-WP/FX-WP, the breeze break is not secured. Therefore, handle the balance with care when transporting, etc.


Models : FZ / FX, FZ-WP / FX-WP

2. Unpacking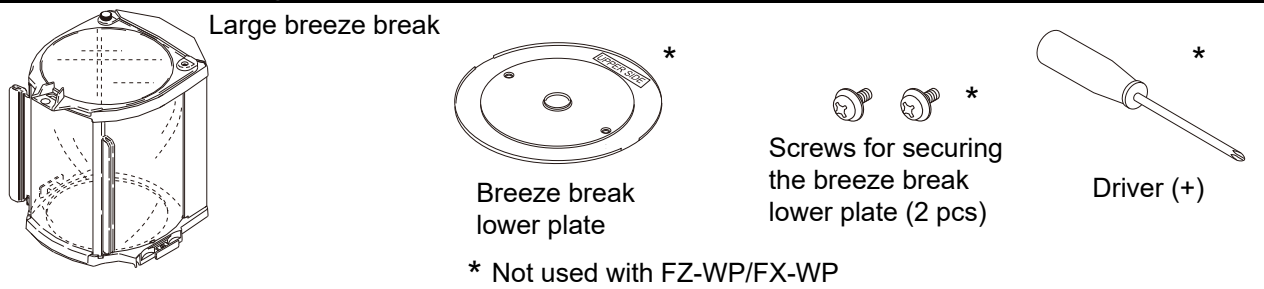


3. Installing the large breeze break

**Note:** For the FZ-WP/FX-WP, steps 1 to 3 do not need to be performed; simply place the breeze break on top.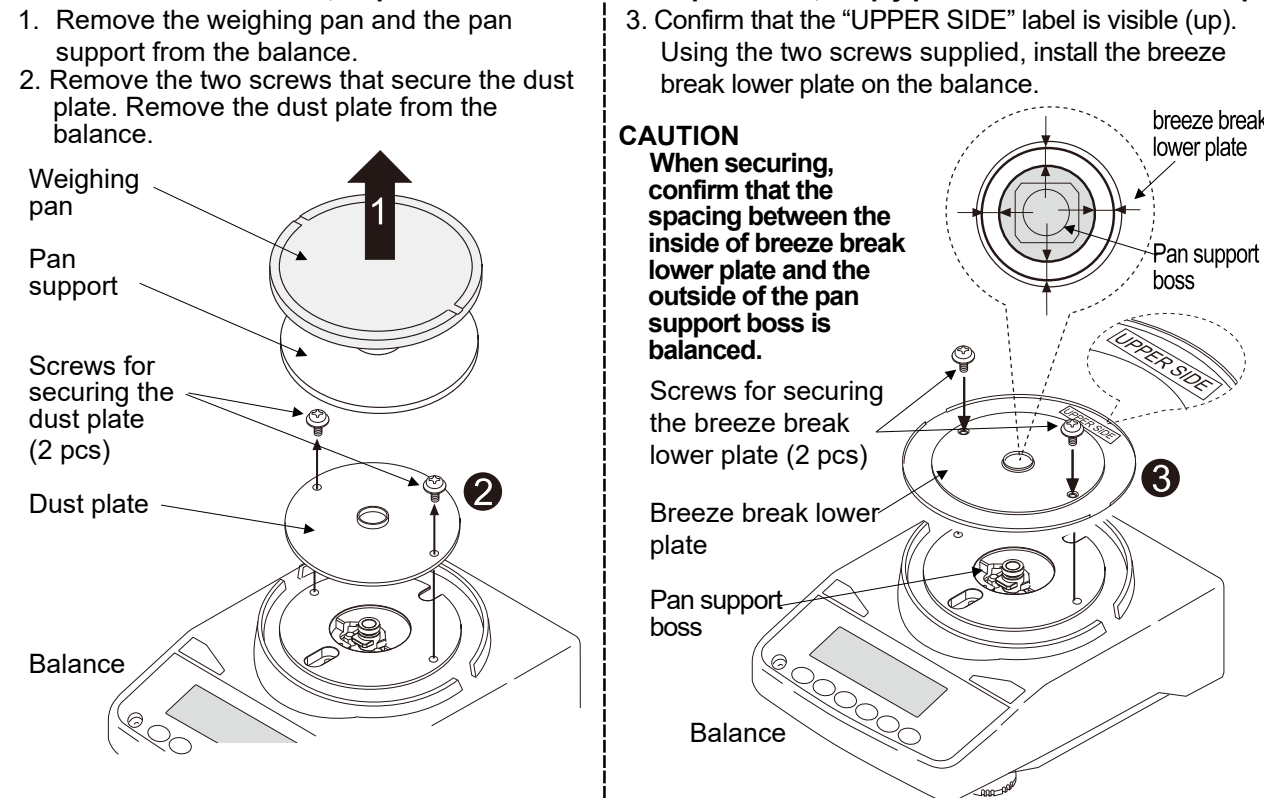


**CAUTION**

- When installing, use care not to apply excessive load to the sensor mechanism (circumference of the pan support boss) of the balance.
- When securing the breeze break lower plate, use the accessory screws provided with the FX-11. (Keep the original dust plate and screws for later reinstallation.)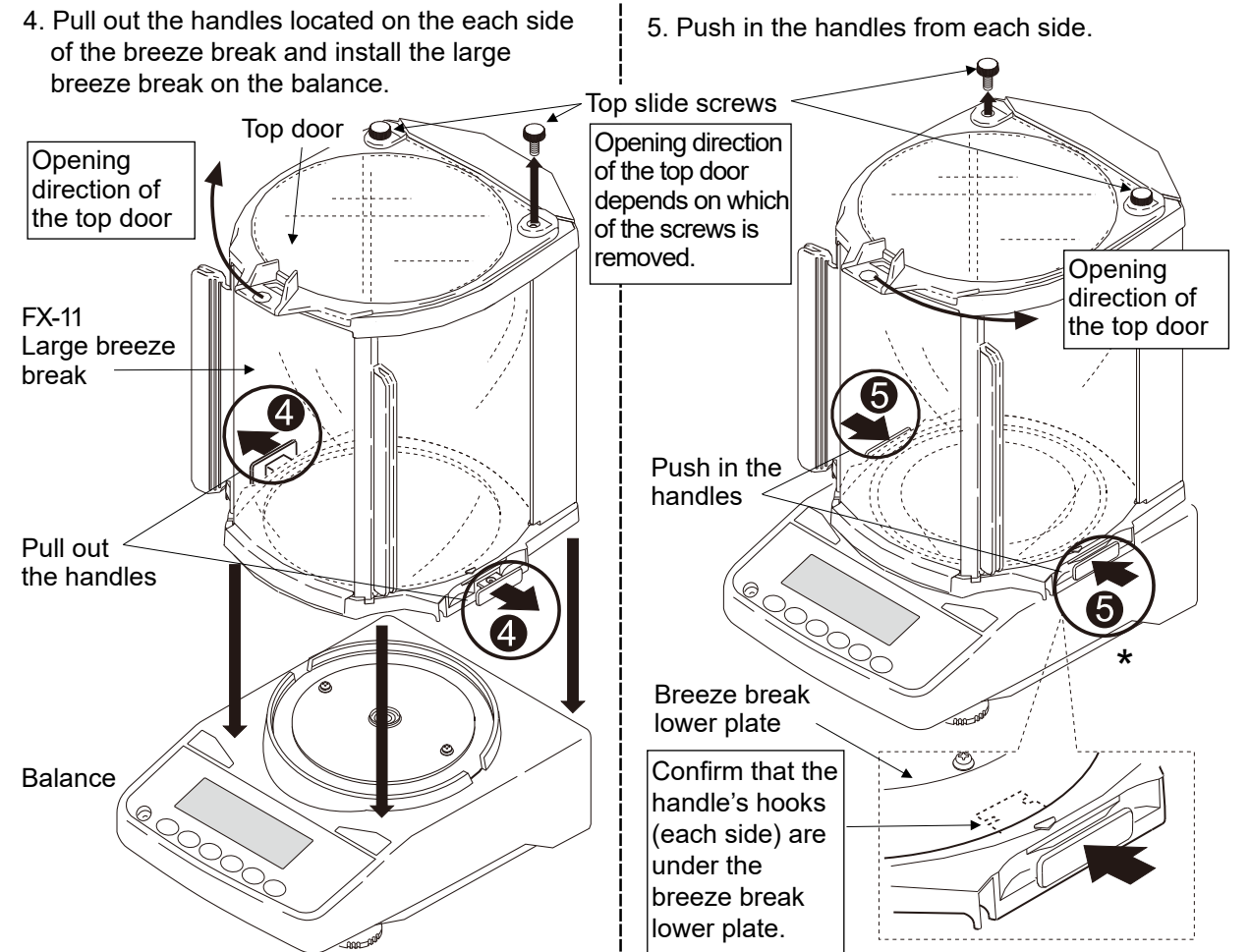

\* When used with FZ-WP / FX-WP the breeze break cannot be secured on the balance whether the handles are pulled out or pushed in.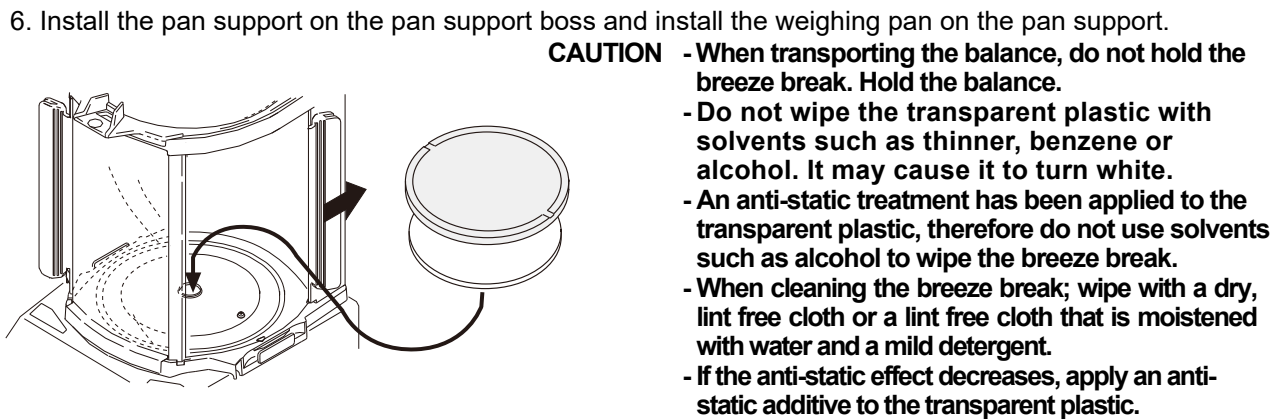

Materials used (Material of the breeze break)		
Transparent plastic: An anti-static treated acrylic	Slide pin: Polyacetal	Other plastic: Antistatic ABS
Screws: steel	Other metal: SUS	